

## Getinge Lancer Ultima labware washing systems

Freestanding series





# Professional cleaning results

## integrated systems,repeatable results

From quality control to general laboratory research, Getinge Lancer Ultima series washers meet space and budget requirements in laboratories performing a range of critical applications where clean laboratories essential.

Typical applications include analytical chemistry, quality control, pharmaceutical and environmental science laboratories, wastewater treatment, food and beverage analysis, public health and consumer product testing facilities.

The exclusive Prolux programmable microprocessor controller commands a full range of prewash, wash, rinse and drying functions through simple touchscreen menus on all Ultima series washers. The convenience of five preset programs for light to heavy loads simplifies operation, while more complex programs can be customized whenever needed to meet specific applications.



Getinge Lancer Ultima model 1300 LX freestanding labware washer, shown with accessory loading rack, sold separately.

## Meeting a variety of laboratory needs

The Getinge Lancer Ultima series freestanding washers include five high-performance models that offer a range of labware loading options using baskets, injector racks or combination racks for the best high capacity processing of multiple labware types in the industry.



910 LX / 1300 LX features	1400 LXP / 1600 LXP features
Prolux control interface	Prolux Plus control interface
Manual fold down door	Manual fold down door
Forced-air drying, injectors, HEPA filtered	Forced-air drying, injectors, HEPA filtered

Capacity comparison					
Model	Volume	Interior (W x H x D)	Positions	Wash levels	Configurations
910 LX	5.7 cu. ft. (160 L)	21.1" x 21.4" x 20.5" (535 x 544 x 520 mm)	3	2	3
1300 LX	6.4 cu. ft. (180 L)	21.1" x 25.4" x 20.5" (535 x 644 x 520 mm)	4	3	5
1400 LXP	10.7 cu. ft. (304 L)	24.6" x 30.4" x 24.8" (626 x 773 x 629 mm)	4	3	6
1600 LXP	14.4 cu. ft. (408 L)	27.7" x 33.8" x 26.7" (703 x 858 x 678 mm)	5	4	12
1600 LXP/HE	17.7 cu. ft. (500 L)	27.7" x 41.3" x 26.7" (703 x 1048 x 678 mm)	5	4	12

Models shown (left to right) 910 LX, 1300 LX, 1400 LXP, 1600 LXP and 1600 LXP/HE.



Getinge Lancer Ultima model 1400 LXP freestanding labware washer, shown with accessory loading rack, sold separately.

Prolux programmable microprocessor controller with intuitive touchscreen user interface.

Control screen shows actual cycle performance, alarm events. Data is captured for documentation and communication.

Preset or user programmed cycles are fully adjustable to permit multiple combinations of prewash, wash, rinse, acid rinse, purified final rinse, HEPA filtered forced air drying, controlled duration, temperature and more.

High efficiency powerful fluid circulation with direct fluid path and rear manifold connections deliver optimal pressure and flow rates.

Chamber walls made in durable AISI 316L stainless steel material.

Injector jet performance provides inside out wash and rinse functions.

Stainless steel top and bottom spray arms and multiple levels of injection washing ensure thorough spray coverage.

Two integrated automatic chemical dosing pumps are standard on all models. Convenient on-board chemical storage beneath chamber saves space.



#### View-In-Process Window

An optional View-In-Process (VIP) window includes an interior LED light for visual verification of the interior load during the cycle.



#### **Cool Surface**

High-performance insulation contains heat and limits energy consumption.



#### **Quiet Operation** Noise reduction

design minimizes sound for quiet operation.

#### Comprehensive collection for any lab

Getinge Lancer Ultima offers a full range of labware washers that address applications at any capacity for highdensity processing of multiple labware types, aiming.to improve throughput. All models are produced to withstand harsh laboratory conditions, minimize water consumption and save energy.

Next to the freestanding washers, the Getinge Lancer Ultima series also includes undercounter washers designed to deliver clean labware faster than other industrial washers. Undercounter washers are available in three configurations and two sizes with generous load capacities.

For more information on undercounter models please refer to the Getinge Lancer Ultima series undercounter brochure.




Getinge Lancer Ultima model 810 LX undercounter washer


5

## Small spaces, impressive capacity

910 LX

Drawings display front and side of unit, with door swing allowance.

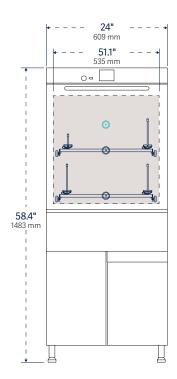

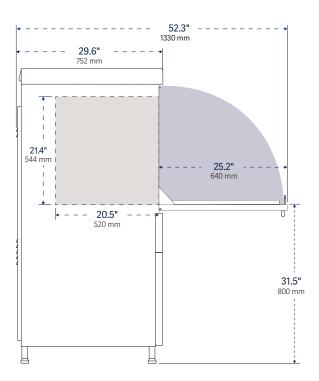

Model	Interior dimensions	Exterior dimensions	Water per fill*	Programs	
	WxHxD	WxHxD	gal (L)	presets	custom
910 LX	21.1" x 21.4" x 20.5" 535 x 544 x 520 mm	24" x 58.4" x 29.6" 609 x 1483 x 752 mm	3.4 (13)	5	35

 $<sup>^{\</sup>star}$  Total water consumption per fill based on reservoir sump and all fluid path fill volumes, care should be taken when evaluating comparative water consumption claims.





Model		Cycle function				Chemical dosing			
	Wash temp. max. °F (°C)	Rinse purified water	Rinse acid	Drying	Automatic pumps	On-Board chemical storage	Level sensors	Water softener	<b>Empty</b> lb (kg)
910 LX	203°F (95°C)		•	Forced-air chamber and injectors; HEPA filtered	2 (max 3)	•	2 (max 3)		315 (143)

■ / 


■ Standard / optional features


7

## Compact footprint, lasting impression

1300 LX

Drawings display front and side of unit, with door swing allowance.

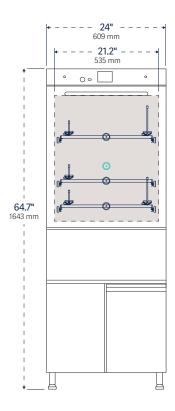


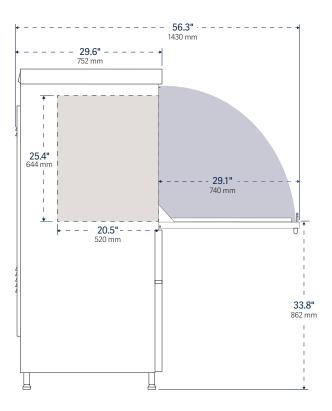

Model	Interior dimensions	Exterior dimensions	Water per fill*	Programs	
	WxHxD	WxHxD	gal (L)	presets	custom
1300 LX	21.1" x 25.4" x 20.5" 535 x 644 x 520 mm	24" x 64.7" x 29.6" 609 x 1643 x 752 mm	4 (15)	5	35

<sup>\*</sup> Total water consumption per fill based on reservoir sump and all fluid path fill volumes, care should be taken when evaluating comparative water consumption claims.



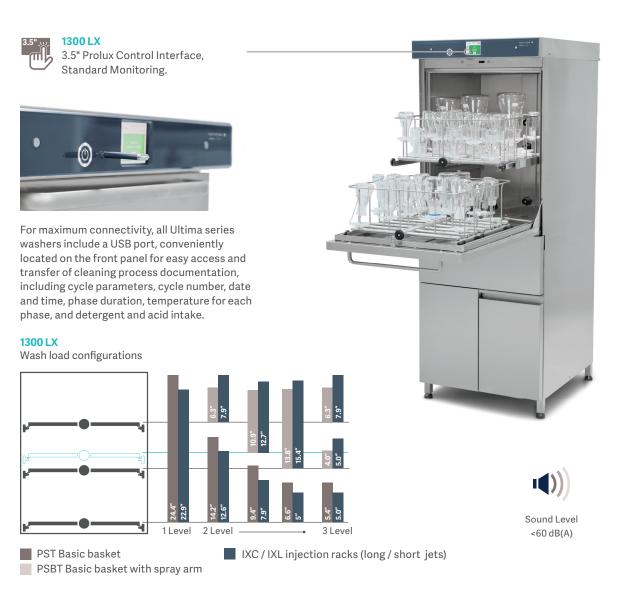

Model		Cycle function				Chemical dosing			
	Wash temp. max. °F (°C)	Rinse purified water	Rinse acid	Drying	Automatic pumps	On-Board chemical storage	Level sensors	Water softener	<b>Empty</b> lb (kg)
1300 LX	203°F (95°C)	•	•	Forced-air chamber and injectors; HEPA filtered	2 (max 3)	•	2 (max 3)		319 (145)


/ Standard / optional features

9

## Space efficiency, multi-level loading

1400 LXP


Drawings display front and side of unit, with door swing allowance.

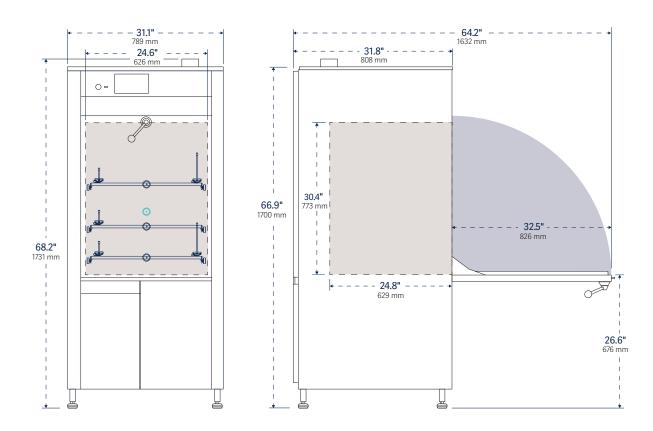

Model	Interior dimensions	Exterior dimensions	Water per fill*	Programs	
	WxHxD	WxHxD	gal (L)	presets	custom
1400 LXP	24.6" x 30.4" x 24.8" 626 x 773 x 629 mm	31.1" x 68.2" x 31.8" 789 x 1731 x 808 mm	5.2-6.6 (20-25)	5	35

<sup>\*</sup> Total water consumption per fill based on reservoir sump and all fluid path fill volumes, care should be taken when evaluating comparative water consumption claims.



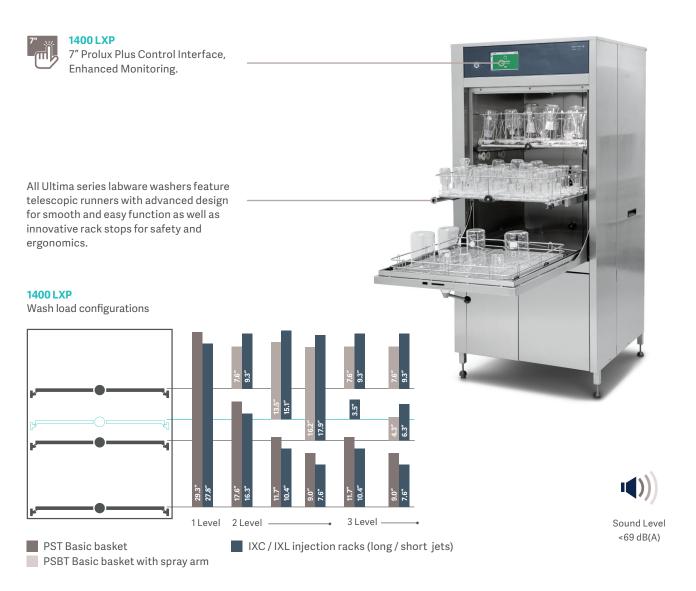

Model		Cycle function				Chemical dosing			
	Wash temp. max. °F (°C)	Rinse purified water	Rinse acid	Drying	Automatic pumps	On-Board chemical storage	Level sensors	Water softener	<b>Empty</b> lb (kg)
1400 LXP	203°F (95°C)	•	•	Forced-air chamber and injectors; HEPA filtered	2 (max 4)	•	2 (max 4)		469 (213)


■ / 
■ Standard / optional features

11

## **Maximum flexibility**

### 1600 LXP

Drawings display front and side of unit, with door swing allowance.

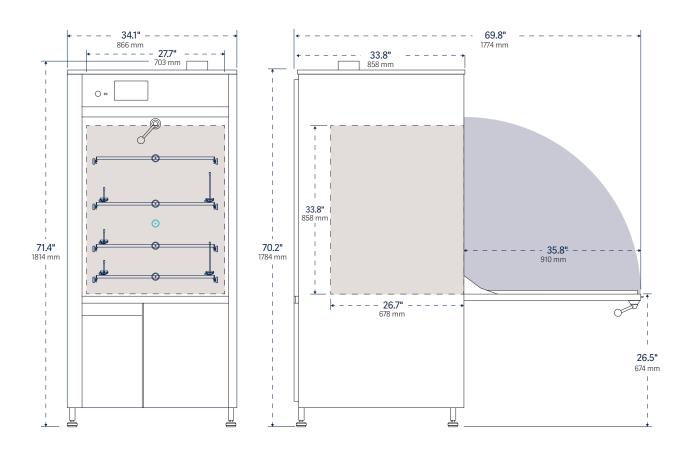

Model	Interior dimensions	Exterior dimensions	Water per fill*	Programs	
	WxHxD	WxHxD	gal (L)	presets	custom
1600 LXP	27.7" x 33.8" x 26.7" 703 x 858 x 678 mm	34.1" x 71.4" x 33.8" 866 x 1814 x 858 mm	7.9-9.2 (30-35)	5	35

 $<sup>^{\</sup>star}$  Total water consumption per fill based on reservoir sump and all fluid path fill volumes, care should be taken when evaluating comparative water consumption claims.





#### 1600 LXP


7" Prolux Plus Control Interface, Enhanced Monitoring.





1600 LXP

Wash load configurations.

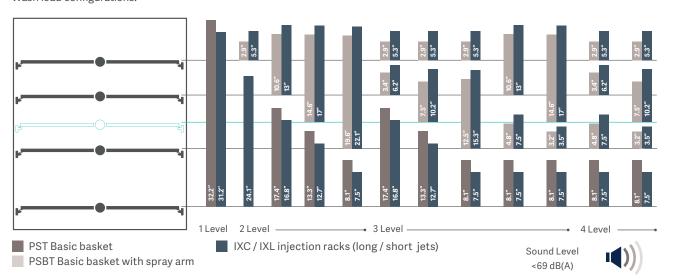

Model		Cycle function				Chemical dosing			
	Wash temp. max. °F (°C)	Rinse purified water	Rinse acid	Drying	Automatic pumps	On-Board chemical storage	Level sensors	Water softener	<b>Empty</b> lb (kg)
1600 LXP	203°F (95°C)		•	Forced-air chamber and injectors; HEPA filtered	2 (max 4)	•	2 (max 4)		474 (215)

■ / 


■ Standard / optional features


13

## **High throughput**

### 1600 LXP/HE

Drawings display front and side of unit, with door swing allowance.

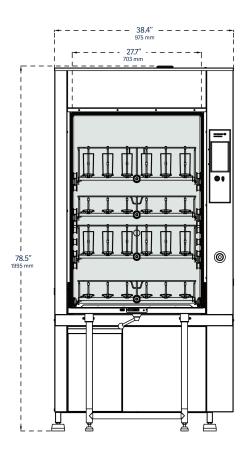

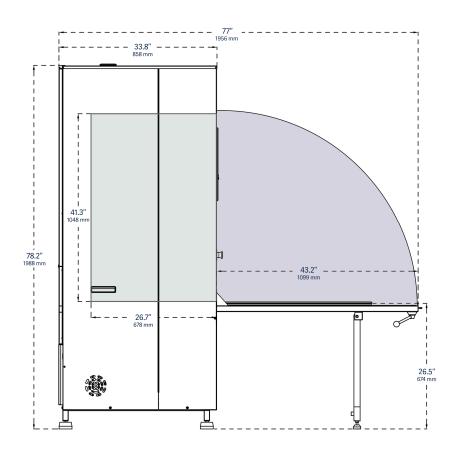

Model	Interior dimensions	Exterior dimensions	Water per fill*	Programs	
	WxHxD	WxHxD	gal (L)	presets	custom
1600 LXP/HE	27.7" x 41.3" x 26.7" 703 x 1048 x 678 mm	38.4" x 78.5" x 33.8" 975 x 1995 x 858 mm	7.9-9.2 (30-35)	5	35

 $<sup>^{\</sup>star}$  Total water consumption per fill based on reservoir sump and all fluid path fill volumes, care should be taken when evaluating comparative water consumption claims.



15



#### 1600 LXP/HE

High throughput chamber with a total loading capacity of 17,7 cu.ft (500 liters)


Include door jacks for safe loading and unloading



#### 1600 LXP/HE

Wash load configurations.

GETINGE LANCER ULTIMA FREESTANDING SERIES

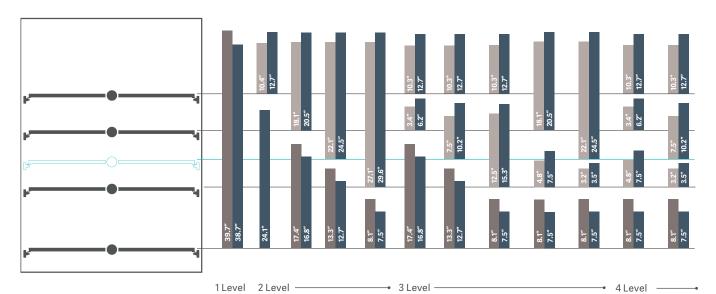

PST Basic basket

IXC / IXL injection racks (long / short jets)

Sound Level

FSBT Basic basket with spray arm

Model	Cycle function				Chemical dosing				Net weight
	Wash temp. max. °F (°C)	Rinse purified water	Rinse acid	Drying	Automatic pumps	On-Board chemical storage	Level sensors	Water softener	<b>Empty</b> lb (kg)
1600 LXP/HE	203°F (95°C)	•	•	Forced-air chamber and injectors; HEPA filtered	2 (max 4)	•	2 (max 4)		529 (240)

☐/☐ Standard/optional features

### Repeatable results

### - with reliability, safety and high throughput

#### **Ergonomic Loading Configurations.**

Unlike oversized washers with less efficient loading patterns, Getinge Lancer Ultima series freestanding labware washers are designed for high-density loading and the most effective use of available interior space. This efficiency maximizes throughput and reduces capital investments, operating costs and the amount of dedicated lab space needed.

- Telescoping load-bearing rails permit extension of racks for easy loading with or without an optional trolley.
- All racks are interchangeable between wash levels thereby minimizing the need for additional accessories.
- Racks for Getinge Lancer Ultima series undercounter models and freestanding models 910 LX, 1300 LX are interchangeable.
- The fold-down door creates a platform for proper rack positioning and more comfortable loading and unloading.

#### Cleaning performance and safety.

Operating systems are designed for performance, efficiency and high throughput.

- An automatic door lock prevents door opening during active cycles and protects the user from heat and chemical exposure.
- Top and bottom stainless steel spray arms are pressurized to provide exterior coverage to the labware.
- Injector rack manifold connections are located at the rear for proper alignment and the most efficient pressure and flow.
- Washing temperature can be set to a maximum of 203°F (95°C) if required for more difficult soils.
- Automatic Peristaltic pumps are self-contained for accurate and repeatable dosing of detergents or acids required by each cycle.
- Multiple DI rinses can be programmed to eliminate interior and exterior spotting.




The front-loading door folds down to create an ergonomically stable platform for proper rack positioning and easier loading, eliminating the need for a trolley to load or unload glassware.

#### Alarms and monitoring

- The high resolution touchscreen color display is easy to read from a distance.
- The intuitive graphic user interface creates a series of visual performance indicators such as active program, time remaining, warning, alarm and cycle completed functions.
- Five preset programs permit one touch operation.
   Additional cycles are programmed from the graphic user interface.
- An integrated USB port allows easy front panel access for process tracking and data logging. Optional process record printouts available through USB or Ethernet ports are managed by the Prolux controller.

Hydraulic design

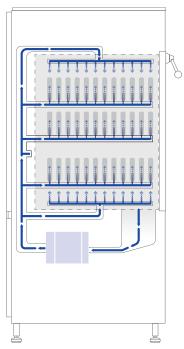

The Getinge Lancer Ultima hydraulic design optimizes pressure and flow for efficient performance at all wash levels.

The combination of a highperformance recirculation pump, straight-line fluid path and center-mounted rack interface assures maximum water flow through spray arms and injectors.

Washer pump is protected from debris by a stainless steel sump screen. The screen lifts out without tools for easy cleaning.

Model 1400 LXP, waterflow.

#### Construction

The sump, spray arms, interior chamber walls and door are fabricated from 316L stainless steel to resist aggressive chemicals and residues.

- Exterior 304L stainless steel panels resist corrosion, simplify cleaning and protect electrical and plumbing components.
- Double-wall insulated assembly confines heat loss and minimizes noise.
- Easy access stainless steel chamber filters extend washer life by protecting pumps and interior plumbing from debris.
- A compact liquid sump and direct fluid path minimizes water volume per fill, lowers overall consumption per cycle and improves sustainability throughout the useful life.

#### **View-In-Process**





An optional dual pane View-In-Process (VIP) window includes an interior LED light for visual verification of the interior load during the cycle. The light is activated by the On/Off switch through the Prolux controller. Tempered safety glass provides excellent sound and heat abatement and enhances operator safety. Factory installed, specify when ordering.

#### Model 910 LX:

13.7" (348 mm) H x 6.8" (173 mm) W **Models 1300 LX / 1400 LXP / 1600 LXP:** 20" (508 mm) H x 6.8" (173 mm) W

\* VIP window not available on model 1600LXP/HE

## Simple, flexible, documented, trusted.

#### Prolux controller.

The all new 3.5" Prolux and 7" Prolux Plus controllers are based on a high-performance PLC microprocessor designed for simplicity, one-touch start, real-time status indicators and intuitive programming options that permit customization over the range of washer operations.

Both the Prolux and Prolux Plus controller integrate a suite of menu screens that support digital functions from cycle selection, process monitoring, warning advisories, audible and visual alarms, system communications and data capture.

#### **Operation**

- Screen menus and graphic user interfaces are designed to simplify setup and operations including active programs, remaining time, warnings, alarms and cycle complete notifications.
- Five preset programs are factory loaded for one-touch control of most routine applications.
- When additional cycle programs are desired, parameters can be changed through secure access to the administration screen where an intuitive graphic user interface displays all options.

#### **Communication and connectivity**

- A front-panel USB port simplifies direct connection to peripheral devices for cycle tracking and data capture.
   The port is covered by a protective hood.
- A real-time process record is printed if optional printer is installed.
- Multiple data ports include Ethernet and RS422 connectors.
- Optional printer documents cycle performance, warnings and alarms.



The Prolux and Prolux Plus (above) controllers include a touchscreen interface with a common icon directory that permits scrolling access to primary screens from the same menu. A set of sub-menu screens display additional icon directories specific to the screen functions.

#### 910 LX / 1300 LX



- 3.5" Prolux color touchscreen interface
- 320 x 240 pixel resolution display
- Five presets
- · Additional custom program options

#### 1400 LXP / 1600 LXP / 1600 LXP/HE



- 7" Prolux Plus color touchscreen interface
- 800 x 480 pixel resolution display
- Five presets
- Additional custom program options



The Prolux control interface features an intuitive menu of primary and secondary screens, configured with icons, to simplify access to all functions. The touchscreen is easily managed with a gloved hand.

#### **Programming**

Programmable cycle parameter selections include:

- Prewash. Up to 3 prewashes, 0 to 30 minutes each, temperatures up to 203°F (95°C), chemical dosing as desired.
- Wash. 0 to 30 minutes, temperatures up to 203°F (95°C), chemical dosing as desired.
- **Acid rinse.** 0 to 30 minutes, chemical dosing, temperatures up to 203°F (95°C).
- Running water rinses. One or more short rinse sequences after the wash or rinse cycle purge the hydraulic circuit of residual detergent and rinsing agents.
- Rinse, DI. Up to 4 rinses, 0 to 30 minutes each. Ambient to 203°F (95°C).
- **Final rinse, DI.** 0 to 30 minutes, temperatures up to 203°F (95°C).
- **Drying.** Up to 90 minutes, adjustable. Temperature, ambient to 230°F (110°C), Prolux Plus models; ambient to 185°F (85°C), Prolux models.
- Cooling. 0 to 30 minutes, adjustable.

#### **Display screens**

Featured display screens are 3.5" screen. Displays for 7" screens are identical.



#### Program Selection.

The Program Selection screen lists the five factory presets, as well as additional presets added through administrator access.



#### Cycle in Process.

A real-time status screen displays time remaining until cycle completion and links to cycle detail screens with the touch of a button.



#### Cycle Completed.

The green Cycle Completed screen confirms completion of the cycle and indicates the washer is safe to unload.



#### Warning Message.

The amber Warning Message screen provides a status report to prompt refill of detergent and rinse agents.



#### Alarm Message.

The red Alarm Message screen confirms the reason for a system shutdown and prompts corrective action needed in advance of a restart.

## **Bioreactor preparation solution**

### - consistent processes deliver repeatable results

The expertise of Getinge and Applikon research and development has established a synergistic process for bioreactor preparation. A solution to prevent cross-contamination, minimize any research downtime and ensure validated outcome through optimal processes and procedures. Getinge offers a specially designed bioreactor wash rack for Getinge Lancer Ultima laboratory washers.

This dedicated wash rack for multi-use bioreactors establishes a highly reproducible system with significant benefits such as:

- Free-up resources to increase laboratory throughput and minimize downtime
- Decreasing operational labor costs compared to manual cleaning processes
- Improving staff safety by removing potential repetitive motions and exposure of harmful substances related to manual procedures.

The relationship between the bioreactor design and laboratory washer assures predictable outcomes.

The defined connection points on the wash rack and bioreactor head plate ensures a clean bioreactor prepared for sterilization.




#### **Baskets**

Small bioreactor assembly parts are placed in the baskets to ensure that these are in place throughout the cleaning cycle.

#### Multiple injector tubes

Reassure best possible cleaning performance for each entry connection on the bioreactor head plate.

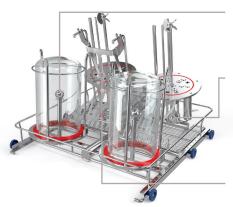

#### **Spray headers**

High flow of process water through the spray headers clean the inside of the bioreactor vessel.

#### **Head plate fixation**

The head plate is fixed in the wash rack. The spray wings within the washer will clean the outside of the head plate.

#### Angled head plate

Minimum water retention on the head plate surface will speed up the drying phase which shorten the overall process cycle.



Getinge Lancer Ultima model 910 LX freestanding labware washer with filled bioreactor wash rack.

#### Bioreactor wash racks for Ultima 910LX and 1300LX

- Cleaning of Getinge Applikon bioreactor size 1-3 liters (0,26-0,79 gal)
- Small bioreactor assembly parts are placed in the baskets to ensure that these are in place throughout the cleaning cvcle.
- Multiple injector tubes reassure best possible cleaning performance for each entry connection on the bioreactor head
- High flow of process water through the spray headers clean the inside of the bioreactor vessel.

#### Bioreactor wash racks for Ultima 1400LXP, 1600LXP and 1600LXP/HE

- Cleaning of Getinge Applikon bioreactor size 1-7 liters (0,26-1,85 gal) but also other branded bioreactors in sizes up to 5
- The bioreactor wash rack includes a base rack to clean various large mouth laboratory glassware and a module (selectable) for the bioreactors docked on the base rack. With this combination we introduce two racks in one.
- Small bioreactor assembly parts are placed in the baskets to ensure that these are in place throughout the cleaning
- Multiple injector tubes reassure best possible cleaning performance for each entry connection on the bioreactor head
- High flow of process water through the spray headers clean the inside of the bioreactor vessel.

#### Configure your bioreactor wash rack, for Ultima 1400 up to 1600LXP/HE models

- 1. Select a base rack, based on your washer model
- 2. Select a module, based on the volume and throughput of your bioreactors
- 3. Place the bioreactors and its assembly parts in the right position








Bioreactors in place



Base rack



Module for 1-7 liter bioreactors

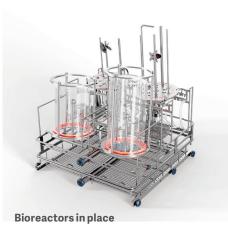

## Adaptable to suit your application

#### Site preparation and utilities

#### **Pre-installation**

Start-up services by factory-authorized technicians include washer set and leveling, pre-installation and final inspections, soft hook-up to utilities (in place by customer), check-out and initial operation. In addition validation documentation and execution options are available; contact us through; sales.lancer@getinge.com

#### Post-installation

Technical support by direct product specialists between 8AM and 7PM Eastern Time EST.

#### **Warranty information**

Getinge Lancer products are supported by a standard industry warranty.

Site preparation	910 LX / 1300 LX / 1400 LXP / 1600 LXP / 1600 LXP/HE
Water supply (hot, cold and DI)	34" hose thread connection, recommended min. flow 5.25 gpm (20 L); lower flow rates may extend cycle times; 29-87 psi (200-600 kPa) water pressure
Drain requirement, vertical standpipe	11/2" (40 mm) diameter, 31" to 35" (800 to 900 mm) above floor.

#### Certifications

UL/IEC 61010-1

#### **Detergents and rinsing agents**

For best cleaning performance there is a selection of detergents and rinse agents formulated to maximize the performance of the Getinge Lancer Ultima labware washer while minimizing water consumption. These cleaning chemicals are made specifically for use in high pressure washers. To order contact sales.lancer@getinge.com.

Getinge LancerClean detergent is available in three solutions formulated for cleaning all different types of contaminants including laboratory soils, "baked on" organic residues, petroleum applications and other hard to clean soils. Getinge LancerClean is ideal for cleaning glassware, surgical instruments and many other labware items.

- LCD-P is a potassium hydroxide based solution that provides excellent cleaning results for all types of contaminants and standard laboratory soils. It is recommended for use in all glassware washers.
- LCD-S is a sodium hydroxide based solution specifically formulated to provide exceptional cleaning for the most difficult types of contaminants.
- LCD-H is a low pH detergent comprised of a sodium silicate solution.

Getinge LCA-A, LCA-C & LCA-P rinse is available in three formulas to meet specific cleaning requirements. Used as a rinsing agent following Getinge LancerClean detergent in the wash cycle, Getinge LCA-A, LCA-C & LCA-P provides excellent results following washing of standard laboratory soils, organics, pharmaceutical based contaminants, petroleum and other difficult materials.

- LCA-A is an acetic acid based neutralizing solution. It is the most commonly used rinsing agent.
- LCA-C is an excellent rinsing agent with a citric acid based solution.
- · LCA-P is a phosphoric acid solution.

#### Some options available

Features		910 LX	1300 LX	1400 LXP	1600 LXP	1600 LXP/HE
Discharge water cooling	Reduces effluent temperature to less than 60°C (140°F) for code compliant transfer to utility drain.					
pH programming	Neutralizes effluent before transfer to utility drain.					
Printer	For hard copy permanent record.					_
Emergency stop	Emergency stop button.					
Sampling valve	Manual sampling valve.					
Conductivity monitoring	Conductivity meter for final rinse verification.					
Pump pressure monitoring	Verification that correct pump pressure is achieved in the water distribution system that secure optimal cleaning performance.					
Watersoftener	Water softener.					
Chemical dosing pump, additional	Additional chemical dosing pump. Up to 1 additional on 910 LX / 1300 LX and up to 2 additional on 1400 LXP / 1600 LXP / 1600 LXP/HE.					
Main Switch	Main power switch, shuts off all power to the washer.	_	_			
Utilities		910 LX	1300 LX	1400 LXP	1600 LXP	1600 LXP/HE
Steam heating	Steam coil to utilize house steam for heating of wash/rinse water.	_	_			
Low pressure kit	Boosts DI water supply input pressure to assure optimum performance.					
Steam condenser	Steam condenser for collapsing exhaust vapor.					
Gravity drain	For quick evacuation of chamber. Requires compressed air and floor drain.	_	_			
Tri-clamp DI inlet	Direct DI water inlet with tri-clamp connection and dry contact for control of customer DI loop valve.	_	_			
Construction		910 LX	1300 LX	1400 LXP	1600 LXP	1600 LXP/HE
Movingrunners	Customized location of rack manifolds to accommodate specific loads.	_	_			
Extended diptube kit	Extended diptubes for use with 33-55 gal (125-208 L) chemical containers.	_	_			
Dry contact	Dry contacts for external communication.					
Seismic restraints	Seismic tie down brackets.					
Stainless steel plumbing	Complete stainless steel hydraulic circuit including main circulation pump	_	_			
View-In-Process window	Includes interior light, permitting visual verification during wash cycles.					_

/ Standard / optional features


 $<sup>\</sup>ensuremath{^{\star}}$  See product specifications for details and additional options to select.

### Common accessories

**Choose from up to 60 different injector racks**, pipette racks and baskets. All racks are interchangeable between the washer lower and upper levels.

- A wide variety of standard inventory baskets and injector racks are available to meet most needs.
- Custom racks and baskets can be fabricated to meet specific needs or special applications.

Please note. All dimensions in the accessories section are in millimeters, unless otherwise stated.

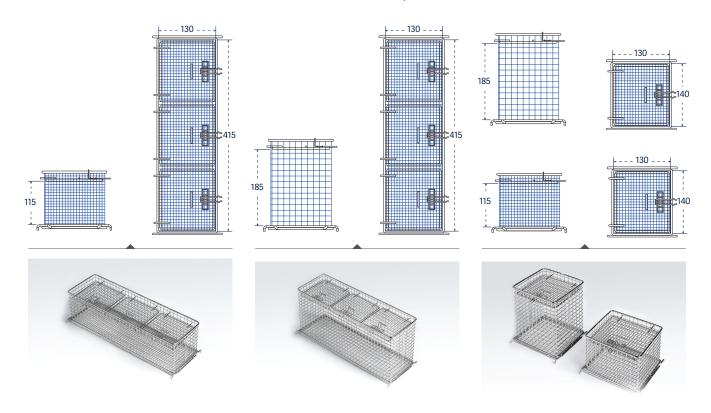

Small mesh basket with cover, long: Catalog No. STCXL

Small mesh basket, tall with cover, long: Catalog No. LTCXL

Small mesh baskets, tall and short: Catalog No. LTC and STC

Small mesh baskets For bottle caps, ferrules, plugs, stoppers and test tubes.	910 LX / 1300 LX / 1400 LXP / 1600 LXP / 1600 LXP/HE	Part number
Mesh basket, short, with cover, 130x118x98 mm	STC	02010037
Mesh basket, short, with cover, 405x118x98 mm	STCXL	02010038
Mesh basket, tall, with cover, 130x118x170 mm	LTC	02010029
Mesh basket, tall, with cover, 405x118x170 mm	LTCXL	02010030





## Accessories

910 LX / 1300 LX





Basic basket Catalog No. PST



Basic basket with spray arm: Catalog No. PSBT



Basket for culture dishes: Catalog No. BP



Support for vials. Catalog VBS




Vials caps support Catalog No. VCS


Baskets Useful for beakers, trays, dishes, dissolution vessels, funnels, instruments and jars.	910 LX / 1300 LX	Part number
Basic basket	PST	AA02600120
Basic basket with spray arm	PSBT	AA02600117


Support holders Useful for culture plates and vials that fit PST and PSBT.	910 LX / 1300 LX	Part number
Support for culture plates to be used with the PST basket	20 PL	AA02250225
Vials bottle support for 2-30 ml vials,	VBS	6007506021
Vials caps support	VCS	6007508131


Basket covers  Accessory basket covers are available to secure small or lightweight labware during the wash cycle.	910 LX / 1300 LX	Part number
Reversible tray lid for PST/PSBT	GC	01030569
Reversible tray lid, half size for PST/PSBT	GC1/2	01030570
Reversible tray, 1/3 of size	GC1/3PS	01030571
Mesh cover for injector rack	MC	01030585
Mesh cover, half size for injector rack	MC1/2PS	01030586
Mesh cover,1/3 of size for injector rack	MC1/3PS	01030587


Short injectors For inside out cleaning of narrow neck glassware, bottles, cylinders, boiling flasks, Erlenmeyer flasks and volumetric flasks.	910 LX / 1300 LX	Part number
Halfrack, 27 short jets, 6 mm diameter	27 IXC 1/2	02600022
36 short jets, 6 mm diameter	36 IXC	02600023
41 short jets, 6 mm BOD rack	41 IXC	02600003
48 short jets, 2 mm diameter	48 IXA	02600026
48 short jets, 6 mm diameter	48 IXC	02600027
64 short jets, 2 mm diameter	64 IXA	02600030
64 short jets, 6 mm diameter	64 IXC	02600032
64 short jets, mixed 2 mm and 6 mm diameter	64 IXA/2	02600031
100 short jets, 2 mm diameter	100 IXA	02600016
100 short jets, 6 mm diameter	100 IXC	02600018
100 short jets, mixed 2 mm and 6 mm diameter	100 IXA/2	02600017

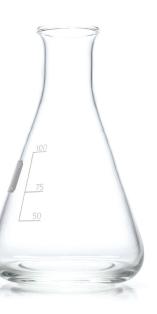


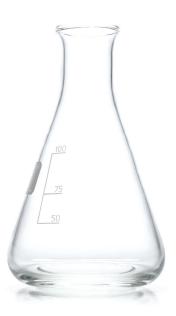


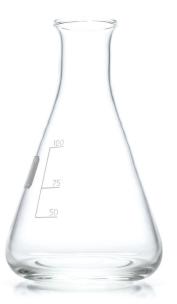


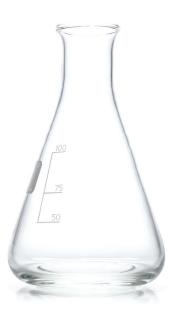

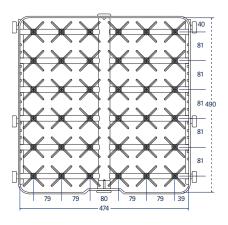


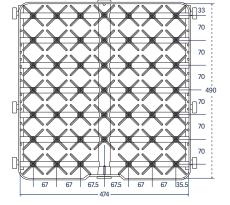

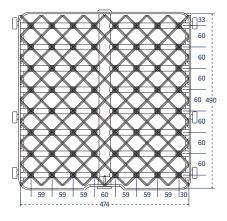

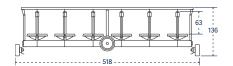


63 136 61.5 136 1.5

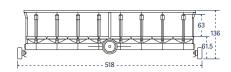

Catalog No. 36 IXC

Catalog No. 48 IXC; IXA

Catalog No. 64 IXA; IXC; IXA/2

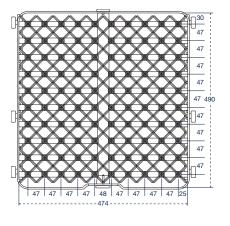



29

IXC: Injector rack, short injectors, 3" (76 mm) tall, jets 6 mm diameter


IXA: Injector rack, short injectors, 3" (76 mm) tall, jets 2 mm diameter


63, 136

Catalog No. 100 IXA; IXC; IXA/2


Long injectors  For inside out cleaning of large narrow neck glassware, bottles, cylinders, boiling flasks, Erlenmeyer flasks and volumetric flasks.	910 LX / 1300 LX	Part number
12 long jets, 8 mm diameter	12 IXL	AA02250224
20 long jets, 8 mm diameter	20 IXL	02600021
36 long jets, 6 mm diameter	36 IXL	02600024
48 long jets, 6 mm diameter	48 IXL	02600028
64 long jets, 6 mm diameter	64 IXL	02600033
100 long jets, 6 mm diameter	100 IXL	02600019

<b>Extra long injectors</b> For inside out cleaning of narrow neck glassware, bottles, cylinders, boiling flasks, Erlenmeyer flasks and large volumetric flasks.	910 LX / 1300 LX	Part number
1 long jet, 12.7 mm diameter, with spray ball	1 IXL	6007506822
2 long jets, 12.7 mm diameter, with spray ball	2 IXL	6007506825
4 long jets, 12.7 mm diameter, with spray ball	4 IXL	6007506828
5 long jets, 12.7 mm diameter, with spray ball	5 IXL	6007506831
9 long jets, 12,7 mm diameter, with spray ball	9 IXL	AA02250223

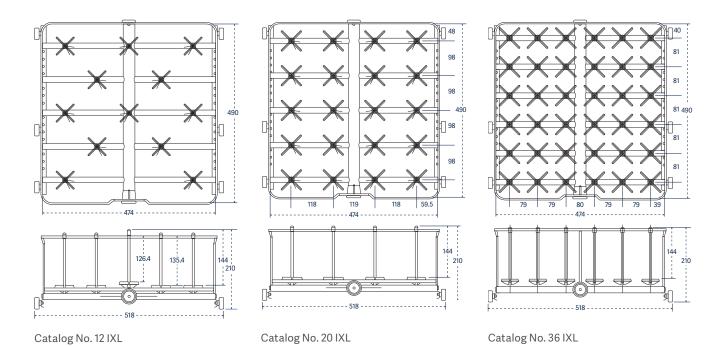


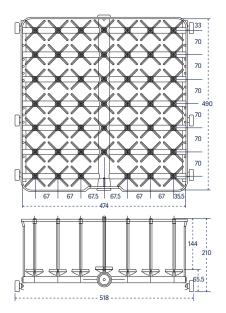

Catalog No. 48 IXL

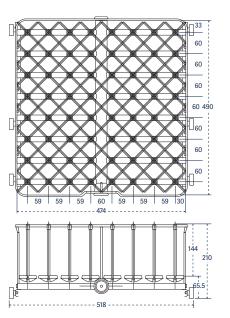

Catalog No. 64 IXL

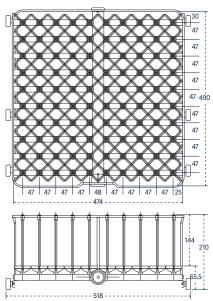

Catalog No. 100 IXL



Catalog No. 1 IXL



Catalog No. 2 IXL



Adjustable clip positions labware on injector for optimal cleaning performance.




Catalog No. 4 IXL



Catalog No. 5 IXL




Catalog No. 9 IXL

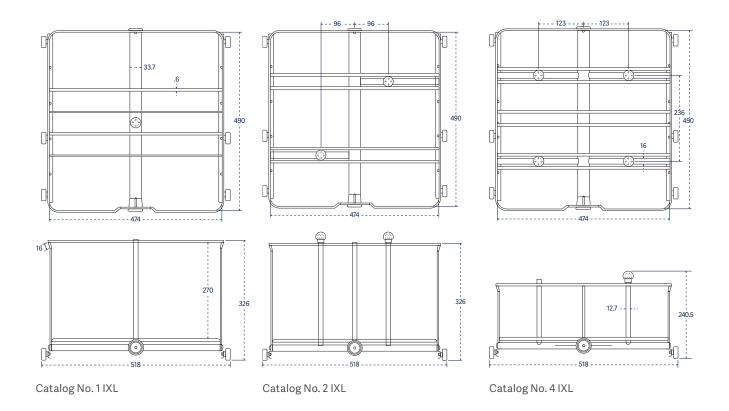

Injector KIT packages	
KIT enable customization of standard injector racks to fit specific needs. Each KIT include 10 complete injectors with their supports and nozzles.	Part number
Long injectors, length of 230 mm and 6 mm diameter	6007500984
Long injectors, length of 171 mm and 6 mm in diameter	6007500890
Short injectors, length of 86 mm and 6 mm in diameter	6007500957
Short injectors, length of 56 mm and 6 mm in diameter	6007501012
Long injectors, length of 171 mm and 8 mm in diameter	6007501011
Needle injectors, length of 86 mm	6007501007



Combination short/long injectors  For inside out cleaning of mixed loads of narrow neck glassware, bottles, cylinders, boiling flasks, Erlenmeyer flasks and volumetric flasks.	910 LX / 1300 LX	Part number
36 short and long jets, 6 mm diameter	36 IXLC	02600025
48 short and long jets, 6 mm diameter	48 IXLC	02600029
48short and long jets, $6mm$ diameter, $2rows2mm$ diameter	48 IXLCA	02600006
64 short and long jets, 6 mm diameter	64 IXLC	02600034
100 short and long jets, 6 mm diameter	100 IXLC	02600020

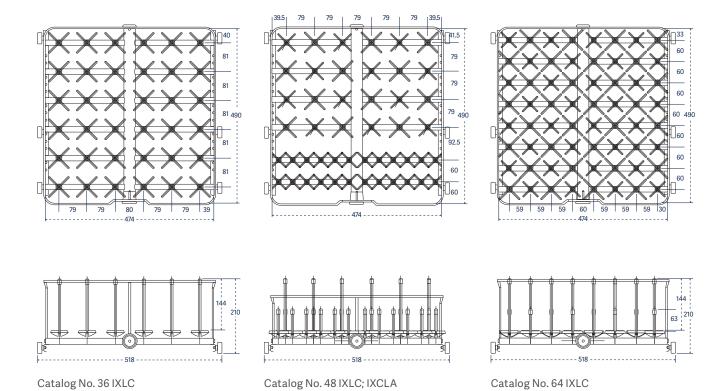

33

<b>Multipurpose racks</b> For a combination of labware and pipettes.	910 LX / 1300 LX	Part number
Basic basket plus 27 short and long jets	PS/27 IXLC	02600039
25 pipettes, angled plus 24 short and long jets	25 PIP/24 IXLC	02600004
50 pipettes, angled plus 34 short and long jets	50 PIP/34 IXLC	02600037
14 short jets plus 4 long injectors angled and perforated	4 EP/14 IXC	02600041
28shortandlongjetsplus4longinjectors, angledandperforated	4 EP/28 IXLC	02600036
One third PS plus 21 short and long jets, plus 25 pipettes	PS/21 IXLC/25 PIP	02600038

Bioreactor rack	910 LX / 1300 LX	Part number
Rack for 1-3 L Getinge Applikon bioreactors	BRWR	AA02600113

Pipette racks		
Accessory pipette racks are configured to accommodate reusable pipettes.	910 LX / 1300 LX	Part number


Grommet style, individually loaded, 100 position

100 PIP









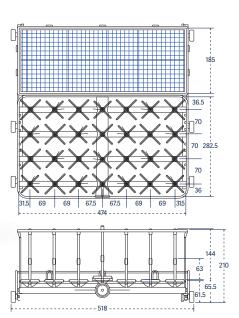

Catalog No. PS/27/IXLC

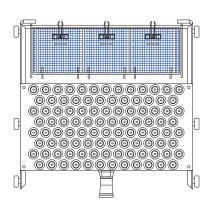

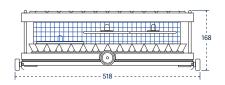

Catalog No. 100 PIP

Storage stand An accessory storage stand offers a convenient place to store accessory racks and baskets when not in use.	910LX/1300LX	Part number
Holds up to 5 racks	RP	01040060

Trolley  The stainless steel loading trolley transfers the load and mates with the washer to simplify loading and unloading. The trolley includes a retainer lip to keep the floor dry from residual water. Includes locking casters.	910 LX	1300 LX
Loading trolley	02530017	02530018



35

GETINGE LANCER ULTIMA SERIES FREESTANDING



## **Accessories**

1400 LXP





Basic basket Catalog No. PST

Baskets

Basic basket

Basic basket with spray arm

Vials bottle support for 2-30 ml vials,

Vials caps support



Basic basket with spray arm: Catalog No. PSBT



Basket for culture dishes: Catalog No. BP

Useful for beakers, trays, dishes, dissolution vessels, funnels, instruments and jars.

Support for culture plates (20 pcs./support max 3 supports) to be



Support for vials. Catalog No. VBS



Part number	
6007504930	

AA02600118

01030417

6007506021

6007508131

1400 LXP	Part number	
20 PL	AA02250225	

VBS

VCS

1400 LXP

PST 14

PSBT 14


20 PL


Basket covers for basic grid baskets Accessory basket covers are available to secure small or lightweight labware during the wash cycle.	1400 LXP	Part number
Reversible tray lid for PST/PSBT	GC 14	02250017
Reversible tray lid, half size for PST/PSBT	GC 1/2 14	02250025

GETINGE LANCER ULTIMA FREESTANDING SERIES 37

Short injectors For inside out cleaning of narrow neck glassware, bottles, cylinders, boiling flasks, Erlenmeyer flasks and volumetric flasks.	1400 LXP	Part number
48 short jets, 6 mm diameter	48 IXC 14	02600059
64 short jets, 6 mm diameter	64 IXC 14	02600063
64 short jets, mixed 2 mm and 6 mm diameter	64 IXA/2 14	02600062
81 short jets, 2 mm diameter	81 IXA	02600067
81 short jets, 6 mm diameter	81 IXC 14	02600068
81 short jets, mixed 2 mm and 6 mm diameter	81 IXA/214	02600066
100 short jets, 2 mm diameter	100 IXA 14	02600054
100 short jets, 6 mm diameter	100 IXC 14	02600053
100 short jets, mixed 2 mm and 6 mm diameter	100 IXA/2 14	02600052



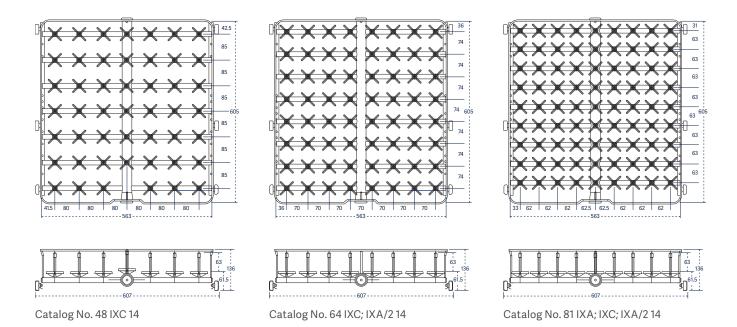

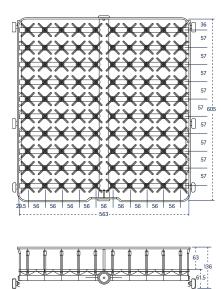


Catalog No. 100 IXA; IXC; IXA/2 14

GETINGE LANCER ULTIMA FREESTANDING SERIES 39

<b>Long injectors</b> For inside out cleaning of large narrow neck glassware, bottles, cylinders, boiling flasks, Erlenmeyer flasks and volumetric flasks.	1400 LXP	Part number
12 long jets, 8 mm diameter	12 IXL 14	02600044
24 long jets, 8 mm diameter	24 IXL 14	AA02600116
36 long jets, 6 mm diameter	36 IXL 14	02600057
48 long jets, 6 mm diameter	48 IXL 14	02600060
64 long jets, 6 mm diameter	64 IXL 14	02600064

Extra long injectors, 305 mm (12") tall  For inside out cleaning of narrow neck glassware, bottles, cylinders, boiling flasks, Erlenmeyer flasks and large volumetric flasks.	1400 LXP	Part number
1 long jet, 12.7 mm diameter, with spray ball	1 IXL 14	6007506823
2 long jets, 12.7 mm diameter, with spray ball	2 IXL 14	6007506826
4 long jets, 12.7 mm diameter, with spray ball	4 IXL 14	6007506829
5 long jets, 12.7 mm diameter, with spray balls	5 IXL 14	6007506832
9 long jets, 12.7 mm diameter, with spray balls	9 IXL 14 DIFF	02600070



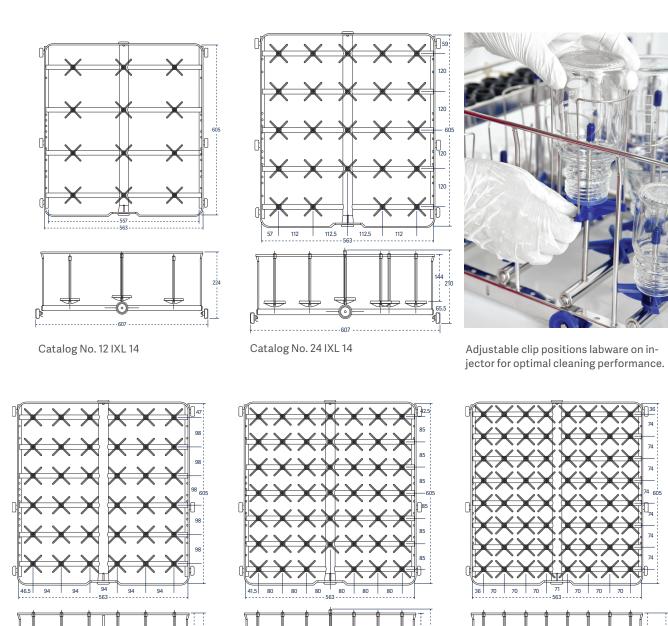

Catalog No. 36 IXL 14 Catalog No. 48 IXL 14 Catalog No. 64 IXL 14

41

Injector rack covers Injector rack covers are available to secure small or lightweight labware during the wash cycle.	1400 LXP	Part number
Injector rack cover full-size	GCI 14	02250023
Injector rack cover half-size	GCI 1/2 14	02250024
Adjustable cross grid	GA 14	02250051









Catalog No. 1 IXL 14

Catalog No. 2 IXL 14


Catalog No. 4 IXL 14

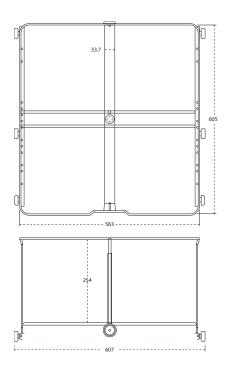

Catalog No. 1 IXL 14

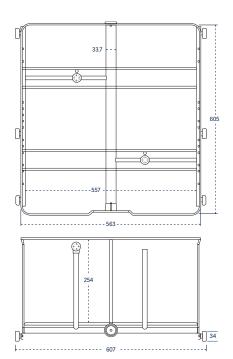



Catalog No. 2 IXL 14

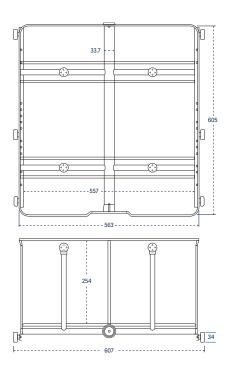


Catalog No. 4 IXL 14


Combination short/long injectors  For inside out cleaning of mixed loads of narrow neck glassware, bottles, cylinders, boiling flasks, Erlenmeyer flasks and volumetric flasks.	1400 LXP	Part number
36 short and long jets, 6 mm diameter	36 IXLC 14	02600058
48 short and long jets, 6 mm diameter	48 IXLC 14	02600061
64 short and long jets, 6 mm diameter	64 IXLC 14	02600065
81 short and long jets, 6 mm diameter	81 IXLC 14	02600069

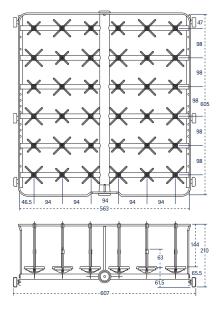

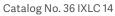

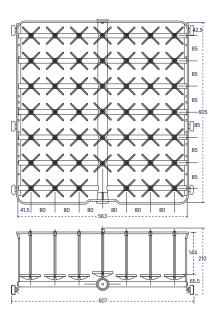

Catalog No. 48 IXL 14

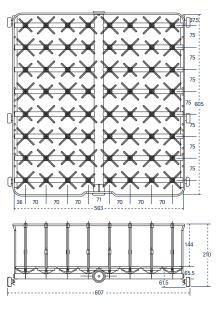



Catalog No. 64 IXLC 14


<b>Multipurpose racks</b> For a combination of labware and pipettes.	1400 LXP	Part number
Basic basket plus 32 short and long jets	PS/32 IXLC 14	02600056
31 pipettes, angled plus 25 short and long jets	31 PIP/24 IXLC 14	02600072
24 short and long jets plus 4 long injectors, angled and perforated	4 EP/24 IXLC 14	02600074


Bioreactor rack	1400 LXP	Part number
Module* for 1-3 L Getinge Applikon bioreactors	BRWRMA	6007506670
Module* for 1-7 L bioreactors	BRWRMB	6007507420
* Basic rack needed for the modules	BRWB 14	6007506671


Pipette racks Accessory pipette racks are configured to accommodate reusable pipettes.	1400 LXP	Part number
Grommet style, individually loaded, 124 position	124 PIP 14	02600012
Pipette rack, canisters (4)	PPI 4B 14	02600075

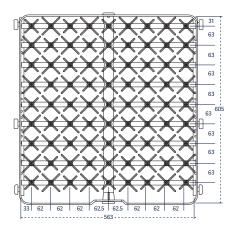


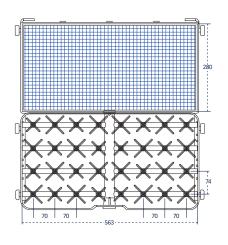

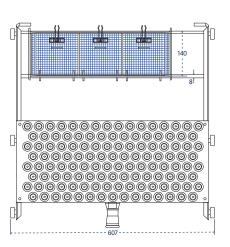

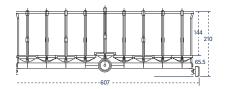

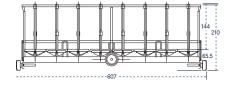

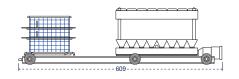

Catalog No. 81 IXLC 14


Catalog No. PS/32 IXLC 14

Catalog No. 124 PIP 14


Injector KIT packages	
KIT enable customization of standard injector racks to fit specific needs. Each KIT include 10 complete injectors with their supports and nozzles.	Part number
Long injectors, length of 230 mm and 6 mm diameter	6007500984
Long injectors, length of 171 mm and 6 mm in diameter	6007500890
Short injectors, length of 86 mm and 6 mm in diameter	6007500957
Short injectors, length of 56 mm and 6 mm in diameter	6007501012
Long injectors, length of 171 mm and 8 mm in diameter	6007501011
Needle injectors, length of 86 mm	6007501007



Storage stand An accessory storage stand offers a convenient place to store accessory racks and baskets when not in use.	1400 LXP	Part number
Holds up to 5 racks	RP 14	02250026
Trolley		
The stainless steel loading trolley transfers the load and mates with the washer to simplify loading and unloading. The trolley includes a retainer lip to keep the floor dry from residual water. Includes locking casters.	1400 LXP	Part number

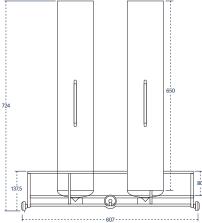



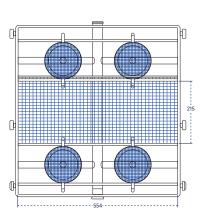



## Accessories

1600 LXP / 1600 LXP/HE





Basic basket Catalog No. PST



Basic basket with spray arm: Catalog No. PSBT



Basket for culture dishes: Catalog No. BP



Support for vials. Catalog No. VBS



Vials caps support Catalog No. VCS


Baskets Useful for beakers, trays, dishes, dissolution vessels, funnels, instruments and jars.	1600 LXP	Part number
Basic basket	PST 16	6007504931
Basic basket with spray arm	PSBT 16	AA02600119
Support for culture plates for PST basket	20 PL	01030417


Support holders Useful for culture plates and vials that fit PST and PSBT.	1600 LXP	Part number
Support for culture plates to be used with the PST basket	20 PL	AA02250225
Vials bottle support for 2-30 ml vials,	VBS	6007506021
Vials caps support	VCS	6007508131


Basket covers Accessory basket covers are available to secure small or lightweight labware during the wash cycle.	1600 LXP	Part number
Reversible tray lid for PST/PSBT	GC 16	01030573
Reversible tray lid for injector racks	GCI 16	01030589
Mesh cover for injector rack	MC	01030585
Mesh cover, half size for injector rack	MC1/2PS	01030586
Mesh cover,1/3 of size for injector rack	MC1/3PS	01030587

Short injectors For inside out cleaning of narrow neck glassware, bottles, cylinders, boiling flasks, Erlenmeyer flasks and volumetric flasks.	1600 LXP	Part number
36 short jets, 6 mm diameter	36 IXC 16	02600096
64 short jets, 6 mm diameter	64 IXC 16	02600100
100 short jets, 2 mm diameter	100 IXA 16	02600088
100 short jets, 6 mm diameter	100 IXC 16	02600090
100 short jets, mixed 2 mm and 6 mm diameter	100 IXA/2 16	02600089
144 short jets, 2 mm diameter	144 IXA 16	02600093
144 short jets, 6 mm diameter	144 IXC 16	02600095
144 short jets, mixed 2 mm and 6 mm diameter	144 IXA/2 16	02600094



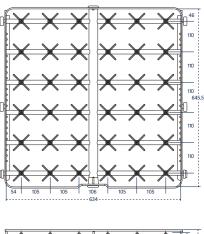


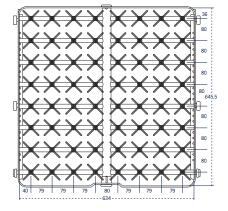


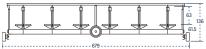


Catalog No. 36 IXC 16

Catalog No.64 IXC 16

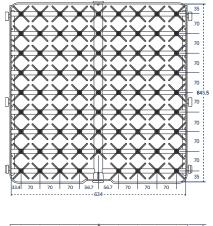

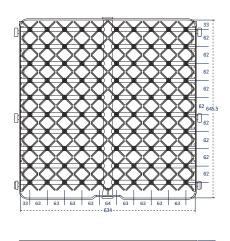

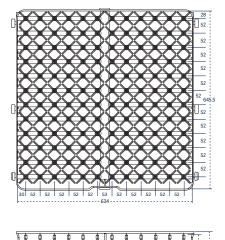


Catalog No. 80 IXC 16


Catalog No. 100 IXA; IXC; IXA/2 16

Catalog No. 144 IXA; IXC; IXA/2 16


<b>Long injectors</b> For inside out cleaning of large narrow neck glassware, bottles, cylinders, boiling flasks, Erlenmeyer flasks and volumetric flasks.	1600 LXP	Part number
12 long jets, 6 mm diameter	12 IXL 16	02600077
36 long jets, 8 mm diameter	36 IXLC 16 D8	02600097
64 long jets, 6 mm diameter	64 IXL 16	02600101
80 long jets, 6 mm diameter	80 IXL 16	02600104
100 long jets, 6 mm diameter	100 IXL 16	02600091

Extra long injectors, 305 mm (12") tall  For inside out cleaning of narrow neck glassware, bottles, cylinders, boiling flasks, Erlenmeyer flasks and large volumetric flasks.	1600 LXP	Part number
1 long jet, 12.7 mm diameter, with spray ball	1 IXL 16	6007506824
2 long jets, 12.7 mm diameter, with spray ball	2 IXL 16	6007506827
4 long jets, 12.7 mm diameter, with spray ball	4 IXL 16 DIFF	6007506830
5 long jets, 12.7 mm diameter, with spray ball	5 IXL 16	02600080
9 long jets, 12.7 mm diameter, with spray balls	9 IXL 16 DIFF	AA02600112

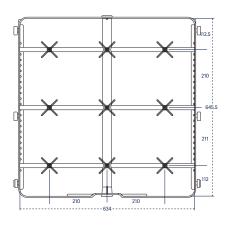

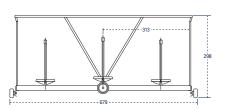


79.5 158 79 158 79.5

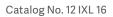





Adjustable clip positions labware on injector for optimal cleaning performance.


Catalog No. 9 IXL 16

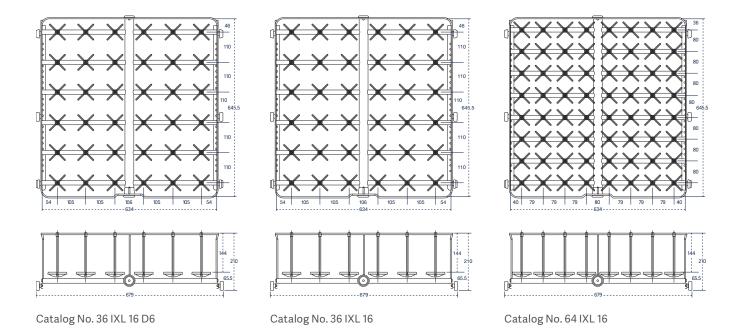


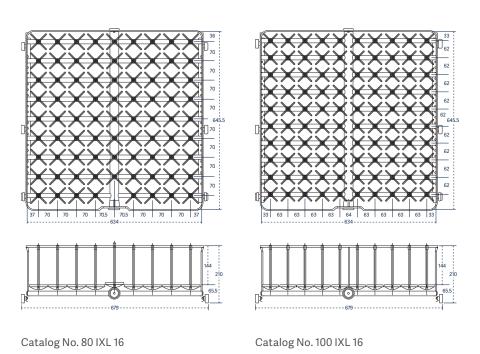


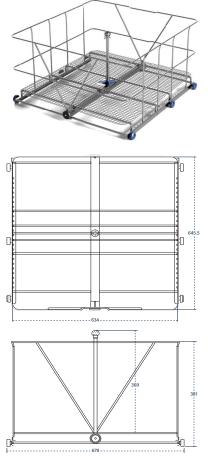


Catalog No. 1 IXL 16

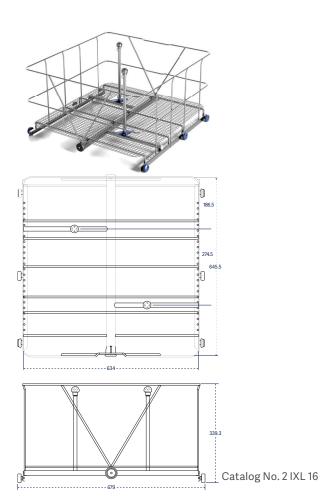


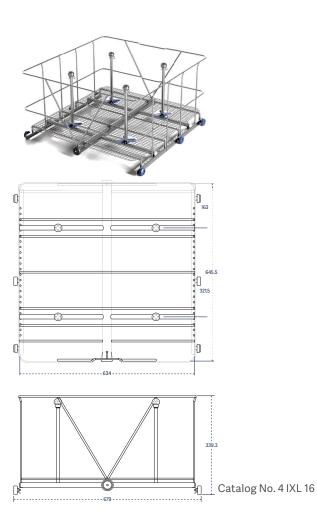


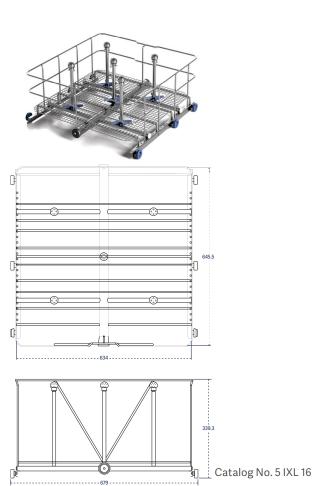



Combination short/long injectors  For inside out cleaning of mixed loads of narrow neck glassware, bottles, cylinders, boiling flasks, Erlenmeyer flasks and volumetric flasks.	1600 LXP	Part number
36 short and long jets, 6 mm diameter	36 IXLC 16	02600099
64 short and long jets, 6 mm diameter	64 IXLC 16	02600102
80 short and long jets, 6 mm diameter	80 IXLC 16	02600105
100 short and long jets, 6 mm diameter	100 IXLC 16	02600092

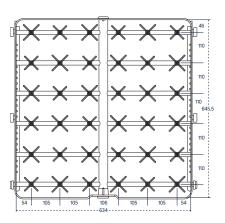


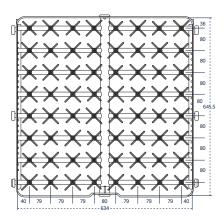

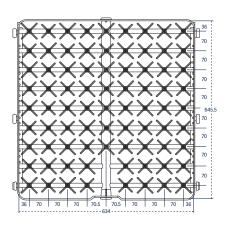

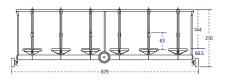

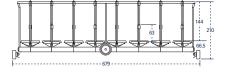

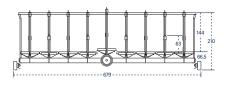

Catalog No. 36 IXLC 16


Catalog No. 64 IXLC 16

Catalog No. 80 IXLC 16

Pipette racks Accessory pipette racks are configured to accommodate reusable pipettes.	1600 LXP	Part number
Grommet style, individually loaded, 105 position	105 PIP 14	02600106

Bioreactor rack	1600 LXP	Part number
Module* for 1-3 L Getinge Applikon bioreactors	BRWRMA	6007506670
Module* for 1-7 L bioreactors	BRWRMB	6007507420
* Basic rack needed for the modules	BRWB 16	6007509429

Storage stand An accessory storage stand offers a convenient place to store accessory racks and baskets when not in use.	1600 LXP	Part number
Holds up to 5 racks	RP 16	01040063

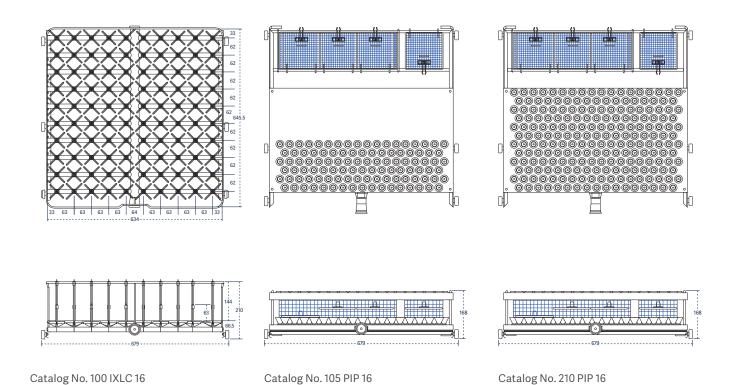

Injector KIT packages	
KIT enable customization of standard injector racks to fit specific needs. Each KIT include 10 complete injectors with their supports and nozzles.	Part number
Long injectors, length of 230 mm and 6 mm diameter	6007500984
Long injectors, length of 171 mm and 6 mm in diameter	6007500890
Short injectors, length of 86 mm and 6 mm in diameter	6007500957
Short injectors, length of 56 mm and 6 mm in diameter	6007501012
Long injectors, length of 171 mm and 8 mm in diameter	6007501011
Needle injectors, length of 86 mm	6007501007



Trolley The stainless steel loading trolley transfers the load and mates with the washer to simplify loading and unloading. The trolley includes a retainer lip to keep the floor dry from residual water. Includes locking casters.	1600 LXP	Part number
Trolley, 4 levels	TN 16	AA02530020



GETINGE LANCER ULTIMA SERIES FREESTANDING 55



Getinge is a global provider of innovative solutions for Life Science companies and institutions, operating rooms, intensive care units and sterilization departments. Based on our firsthand experience and close partnerships with Life Science companies, clinical experts, healthcare professionals and medtech specialists, we are improving everyday life for people – today and tomorrow.

This document is intended to provide information to a North American audience.

Getinge USA Sales · LLC 1150 Emma Oaks Trail Suite 140 · Lake Mary FL 32746 · USA Phone: 1-800-332-1855 · E-mail: Sales.lancer@getinge.com